

Rodent and Small Animal Operating Tables



Heated Rodent Operating Table

- Available as stand alone table or as complete homeothermic heated table
- Uniformly heated with temperature regulation

The Heated Rodent Operating Table is a sturdy piece of equipment that will stand up to daily use. The entire top portion of the table is mounted on a heavy base by means of a

sturdy ball and socket joint. The socket allows the table top to be swiveled and tilted into any convenient position and then locked securely in place.

The stand alone table has a warming chamber under the table top and contains a small, metal-sheathed 40 watt heater. An energy regulator controls the table top surface varying the heating cycle when necessary. The regulator is calibrated from "full" to "off" through ten marked points. Uniform heat distribution is maintained over the entire surface by an insulation barrier located between the woven resistance mat and stainless steel top.

The Homeothermic Rodent Operating Table includes a rectal thermometer-sensing probe and a sensitive control unit. The system continuously monitors the animal's core temperature during surgery and, through the controller, automatically adjusts the table temperature. For complete details on the probes and control unit, see page D2.

These operating tables are supplied with a cork mat that has a cutout in the center for holding the animal. Four stainless steel 6.5 mm (1/4 in) diameter rods and four tying cleats are supplied with the Table. Each 20.4 cm (8-1/16 in) rod is mounted on a simple slotted clamp with a locking screw. The table is mounted on a heavy base fitted with non-slip feet. The top measures 20 x 26 cm (8 x 10 in). It weighs 4.3 kg (9-1/2 lb).

Order #	Product
PY2 50-1205	Heated Rodent Operating Table, 115 VAC, 60 Hz
PY2 50-1213	Heated Rodent Operating Table, 230 VAC, 50 Hz
PY2 50-1221	Replacement Cork Mat
PY2 50-7226-F	Homeothermic Rodent Operating Table with Flexible Probe, 115 VAC, 60 Hz
PY2 50-7227-F	Homeothermic Rodent Operating Table with Flexible Probe, 230 VAC, 50 Hz
PY2 50-7152	Heated Table Only, for use with Homeothermic Control Units PY2 50-7129 and PY2 50-7137



Portable Down-Draft Small Animal Operating Table

- Captures anesthetic gases
- Removable stainless steel work surface
- Autoclavable
- Specifically designed anesthetic chamber

This Down-Draft Table provides an effective way to decrease operator exposure to waste anesthetic gases during rodent procedures without altering the depth of anesthesia. Optional filters for anesthetics, formalin, organic vapors, hazardous and nuisance dusts help comply with laboratory standards and regulations. Captured air is sent either to the filter located within the unit via a blower fan or to an existing exhaust ventilation system.

The Down-Draft Table includes an anesthetic chamber specially designed for use with this table. Accessories include an anodized aluminum cart with optional drop down shelves, storage cabinet and light fixture. Order Potassium Permanganate Filters for formaldehyde use, Carbon Filters for anesthetics use, and Carbon/Potassium Permanganate 50/50 Filters for both formaldehyde and anesthetics applications.

Dimensions

Down Draft Table, H x W x D	45.7 x 61.0 x 76.2 cm (18 x 24 x 30 in)
Work Surface, W x D	76.2 x 43.2 cm (24 x 17 in)

Order #	Product
PY2 72-6201	Portable Down-Draft Table with Blower
PY2 72-6202	Portable Down-Draft Table with Vent
PY2 72-6203	Accessory Package (Folding Shelves, Cabinet and Light)
PY2 72-6204	Folding Shelves only
PY2 72-6205	Anodized Aluminum Cart
PY2 72-6206	Potassium Permanganate Filter, pkg. of 2
PY2 72-6207	Carbon Filter, pkg. of 2
PY2 72-6208	Carbon/Potassium Permanganate Filter 50/50, pkg. of 2
PY2 72-6209	HEPA Filter, pkg. of 1



Connect With Us

