

## Fluovac Masks



Harvard Apparatus are pleased to introduce the new Fluovac mouse mask; a single arm coaxial mask for the delivery and scavenging of anaesthetic gasses, utilising the Fluovac system.

and obtribute guesses, attribute the ride vas ejecenn		
Order No#	Description	
34-1379	Fluovac mask size 0.5 single arm (mice)	
34-0389	Fluovac mask size 1 dual arm (rat)	
34-0390	Fluovac mask size 2 single arm (G. Pig)	
34-0391	Fluovac mask size 3 single arm (rabbit)	

## Scavenged Anaesthetising Chambers



For use in conjunction with the fluovac. Easy to use –simply plug in a size 1 mask into the socket and turn on the fluovac and anaesthetic gas supply. When the animal(s) are anaesthetised the inner chamber slides forward allowing the gas supply to be diverted and the chamber scavenged. During this process fresh air is drawn into the inner chamber to help scavenge any residual anaesthetic.

Order No #	Description	Dimensions H x W x D
34-0455	Scavenged Anaesthetising chamber large	110 x 160 x 260mm
34-0456	Scavenged Anaesthetising Chamber Small	98 x 106 x 130mm

## Fluovac Anaesthetising Chambers



Available as chamber only, or as full gas scavenging systems, including all necessary tubing, connectors and valves. The complete fluovac anaesthetising chamber utilises two directavalves; one is placed on the outlet tube mount of the vaporiser so that anaesthetic gas can be directed to either a face mask or induction box, and the other is placed on the fluovac so that gas can be scavenged from mask or chamber as required.

Order No #	Description	Dimensions H x W x D
34-0460	Anaesthetising Chamber Only Large	200 x 200 x 305mm
34-0462	Anaesthetising Chamber Only Small	134 x 128 x 170mm
34-0458	Large chamber gas scavenging system	200 x 200 x 305mm
34-0461	Small Chamber gas scavenging system	134 x 128 x 170mm

## Fluovac Tube Restrainer



Used for induction of anaesthesia if the patient cannot be held at the Fluovac mask. Simply plug the Fluovac size 1 mask into the restrainer and insert your animal. The animal is then supplied with anaesthesia and the excess is scavenged. Internal diameter 57mm, length is adjustable up to 200mm.

Order No #	Description
34-0457	Fluovac tube restrainer



Connect With Us











