


Pipette fillers



- made of red rubber (natural caoutchouc)
- for graduated and volumetric pipettes
- control of the functions by squeezing the appropriate valves between thumb and forefinger
- individually packed in foil with short instruction for use
- 100 pieces in a carton (multipack)

| Cat. No |           |                                                                                                                | Unit |
|---------|-----------|----------------------------------------------------------------------------------------------------------------|------|
| 5900000 | Standard  | for pipettes up to approx. 20 ml capacity and 5-8 mm diameter, 3 valves                                        | 1    |
| 5900001 | Universal | for pipettes up to 100 ml capacity and 5-9 mm diameter, 3 valves                                               | 1    |
| 5900002 | Flip (TM) | for pipettes up to 100 ml capacity and 5-8 mm diameter, 2 valves with detachable automatic valve easy cleaning | 1    |

Working instruction Flip (TM)

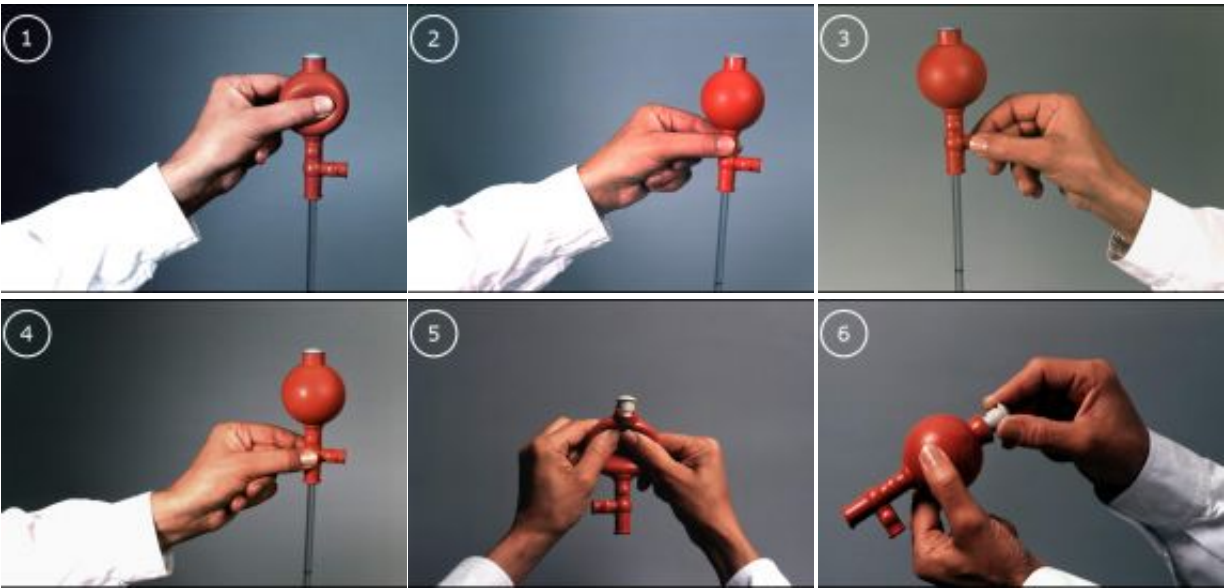

1. Evacuate via the new automatic valve
2. Intake by pressure on operating point
3. Drain by pressure on operating point
4. Blow out the pipette (no valve actuation necessary)
5. Remove the valve for easy inside cleaning
6. Replace the valve



Connect With Us

