


Vacuum Stopping Adapter

Catalog Number: 7593000

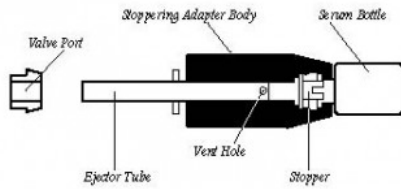

Vacuum Stopping Adapter connects to a valve port for manual stopping of a 20 mm corkage serum bottle under original vacuum. The stopper and serum bottle are inserted into the adapter body and lyophilization begins. After the process is completed, the researcher slides the stopper into position so when the valve is turned from vacuum to vent the serum bottle is tightly sealed by the stopper. The ejector tube is then used to force the serum bottle out of the adapter.

Weight: 1.0 lbs

Weight metric: 0.5 kg

Applicable Products: Serum Bottles