

iWorx RA Recorder

Hardware Manual



iWorx RA Recorder (IX-RA-834)

Overview

The iWorx RA Recorder is the most advanced and feature-rich physiological data acquisition system on the market. It comes standard with eight general purpose input channels, a low voltage and high voltage stimulator, eight digital inputs and outputs, a built-in barometric pressure sensor, and four iWire™ inputs. The iWorx RA exhibits the high resolution, low noise, and flexibility required for a variety of research applications including cardiac function, hemodynamics, isolated tissue and organ studies, exercise physiology research, metabolic function studies, neuroscience and both oocyte and epithelial voltage clamp applications. Incorporating innovative iWire serial interface technology and advanced LabScribe data acquisition and analysis software, the iWorx RA Recorder accommodates a wide range of sensors, transducers, and other devices.

Included Items

- RA Recorder (IX-RA-834)
- USB cable (C-USB)
- Universal 12 VDC power supply (A-UP12-DC-25-1A)
- LabScribe data acquisition software and Quick Start guide (LS-20)

Front Panel



iWorx RA Recorder Front Panel

A1 through A4: BNC connectors for single-ended transducers.



iWorx RA Recorder

A5 through A8: Each channel is equipped with a transducer amplifier to allow connection of virtually any physiologic transducer via a DIN8 connector.

iWire Connectors: Accept up to four serial iWire interfaces including the iWire-B3G, iWire-BIO4, iWire-BIO8, and iWire-ECG12. The iWire-B3G interface can record up to four channels of data. Three of the channels are isolated biopotential amplifiers capable of recording ECGs, EMGs, EOGs, EGGs, and EEGs, while the fourth is a dedicated GSR amplifier (used with the C-ISO-GSR sensor). The iWire-BIO4 and iWire-BIO8 include four or eight biopotential amplifiers respectively. The iWire-ECG12 is capable of recording a 12-lead human electrocardiogram.

S1 and S2: Low voltage stimulators. Parameters for the stimulators, such as pulse width, frequency and amplitude, may be changed on the fly using controls located in the LabScribe software toolbar. Standard protocols include Pulse, Train, Step, Triangle, Ramp, and Custom. Connected via BNC connectors.

HV Stimulator: High voltage isolated stimulator. Suitable for human striated muscle studies. Connected via HV safety connectors.

Rear Panel



iWorx RA Recorder Rear Panel

Power switch

DC Power connector: Accepts 12VDC, 1.5 A DC converter.

Ground: Accepts a banana plug.

EM1 and EM2: Accept the Event Marker (EM-220) and certain other transducers (including the FRS-220 Foot Reaction Switch and the PHRM-100 Polar Heart Rate Monitor receiver).

Sync-Out, Sync-In: To synchronize recordings across multiple RA Recorders so that channels from multiple devices can be sampled simultaneously. BNC connectors.

USB connector

Digital Input/Output 1 and 2: Up to eight digital inputs and outputs are available to monitor and control external TTL devices like pumps and valves.




iWorx RA Recorder

SPECIFICATIONS	
BNC INPUTS (A1-A4)	
Number of inputs	4
Input range	±10 VDC
Resolution	16 bit
Connector	BNC cable
DIN8 TRANSDUCER INPUTS (A5-A8)	
Number of inputs	4
Input range	±10 VDC
Resolution	16 bit
Isolation	No
Excitation	±5 VDC, 100 mA
Connectors	DIN8 (see figure below for pin configuration)
Gain	Programmable with input resistor
HIGH VOLTAGE STIMULATOR OUTPUT	
Connectors	HV Safety
Output range	0-1 mA
Compliance	100 V
Max ON time	10 ms
LOW VOLTAGE STIMULATOR OUTPUTS (S1 and S2)	
Resolution	16 bit
Connectors	BNC
Output range	±15 VDC at 35 mA
Modes	Pulse, Train, Constant, Step, Ramp, Triangle, Custom
DIGITAL INPUTS and OUTPUTS	
Input	8 independent lines, TTL input, 1 Megohm input impedance, 5 V maximum
Output	8 independent lines, TTL output level, 24 mA maximum load per line
A/D CONVERTER	
Sampling speed	200k Hz aggregate
Interface	USB 1.1, 2.0, full speed
GENERAL	
Power	12 VDC, 1.5 A
Enclosure and Dimensions	Plastic; 23 cm W, 15 cm D, 6.5 cm H
Warranty	3 years
Software	LabScribe recording and analysis software



iWorx RA Recorder

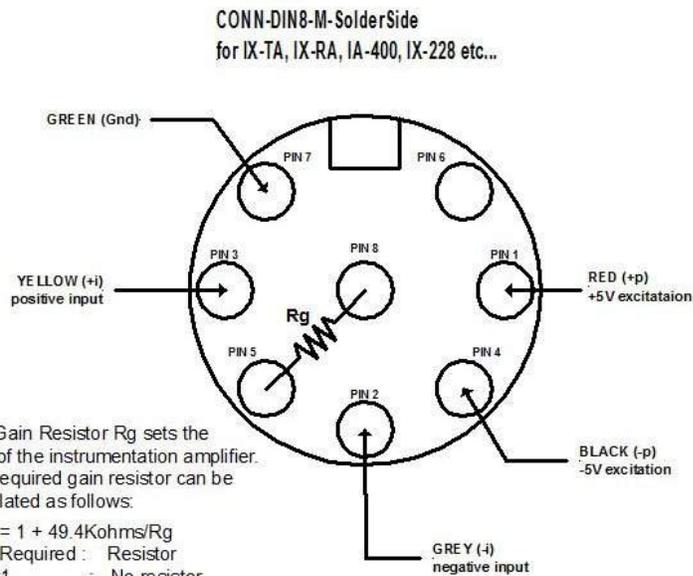

DIN8 transducer connector pin configuration

How to Use the iWorx RA Recorder

The iWorx RA Recorder is controlled by LabScribe recording and analysis software. LabScribe has an intuitive interface for setting up acquisition screens, calibrating signals, and analyzing data on up to 128 channels simultaneously at sampling rates as high as 200K samples/second.

To set up and start to use the iWorx RA Recorder:

- 1) Download LabScribe from www.iworx.com or load the LabScribe software from the provided DVD. Be sure to install the software for the appropriate computer platform (Windows or Mac OS) before connecting the hardware for the first time. In order to download software from the iWorx website, you will need to first register as a user.
- 2) Plug in and switch on the iWorx RA Recorder. If you will be running an experiment that uses an iWire device, connect the device before turning on the RA Recorder.
- 3) Double-click on the LabScribe shortcut to open the software. You should see a message indicating that the iWorx RA Recorder has been recognized by the software.
- 4) There are a number of advanced software modules available for LabScribe. These include modules for research involving intravascular blood pressure, ventricular pressure-volume loops, metabolic function, and ECG analysis.
- 5) Connect the desired transducers to the appropriate channels.



iWorx RA Recorder

- 6) If you need to connect an iWire device once the RA Recorder is already running, it must be turned off, the device connected, and then the RA Recorder turned back on.
- 7) Channels can be configured manually by using the LabScribe Preferences, accessed in the Windows Edit menu, and the Macintosh LabScribe menu. Here channels can be turned on or off, pre-set filters can be selected, and computed channels can be configured. Views can be configured, and stimulator parameters can be set. See the LabScribe User Manual for details.
- 8) Some channel parameters can also be changed from the individual channel menus in the recording window.
- 9) If you are using the Low Voltage Stimulator or the isolated High Voltage Stimulator, connect your stimulating electrodes to the appropriate stimulator and use LabScribe to control the stimulator as instructed in the LabScribe manual.



Connect With Us

